


# B.1.1 Detecting $\gamma$ -Radiation

SG6111

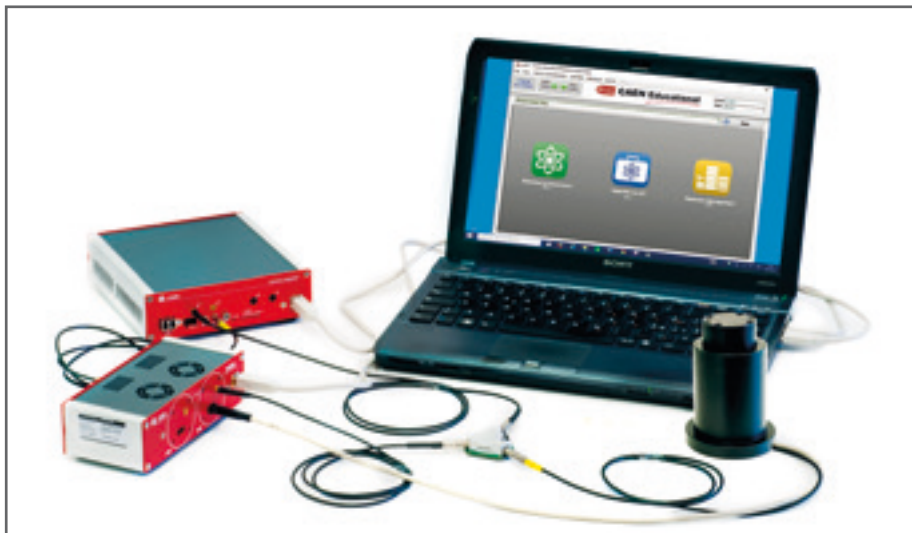

| Dedicated kit                           |     |
|-----------------------------------------|-----|
| Description                             | pp. |
| <b>SP5600C</b><br>Educational Gamma Kit | 179 |



## Requirements

Gamma Radioactive Source 

### Difficulty



### Execution Time



### Data Analysis

NO

### Radioactive Sources

YES

## Equipment

### SP5600C - Educational Gamma Kit

| Model       | SP5600                                                                              | DT5720A                                                                             | A315                                                                                | SP5606                                                                              | SP5607                                                                               |
|-------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|-------------------------------------------------------------------------------------|--------------------------------------------------------------------------------------|
| Description | Power Supply and Amplification Unit                                                 | Desktop Digitizer 250 MS/s                                                          | Splitter                                                                            | Mini-Spectrometer                                                                   | Absorption tool                                                                      |
|             |  |  |  |  |  |
|             | p. 190                                                                              | p. 190                                                                              | p. 192                                                                              | p. 192                                                                              | p. 193                                                                               |

## Purpose of the experiment

*Gamma radioactivity detection by using a system composed of a scintillating crystal coupled to a photon detector.*

See the  
Application



## Fundamentals

Gamma rays interact with matter by three processes: Compton Scattering, Photoelectric Effect and Pair Production (whenever the energy exceeds the 1.022 MeV threshold corresponding to the  $e^+e^-$  rest mass). The cross section of each process depends on the energy of the gamma ray.


The Compton Effect is the inelastic scattering between the incoming photon and an atomic electron. In the Photoelectric Effect, the incident gamma ray transfers all of its energy to a bound electron which acquires a kinetic energy equal to the incoming gamma energy decreased by the binding energy.

These processes convert, totally or partially, the gamma ray energy into kinetic energy of electrons (or positrons, in case of pair production). The interaction of the charged particles with the atomic and molecular systems of the medium results in excited states whose decay, possibly mediated, leads to light in the visible or UV region, eventually detected by the light sensor. A wide range of scintillator products is available today, differing for the light yield, the material properties, the time characteristics of

the scintillation light and, last but not least, cost. The choice of the scintillator is essentially dependent on the specific targeted application.

### Carrying out the experiment


The selected scintillator crystal shall be coupled to the SiPM in the SP5607, through a thin layer of index matching grease to maximize the light collection. In order to avoid saturation, the output of the SiPM is divided using the A315 splitter: one branch is connected to the DT5720A and will be digitized. The other branch will be amplified by the SP5600 module, generating the trigger for the integration signal by the on-board leading edge discriminator or simply counting the pulses induced by the detected gamma ray

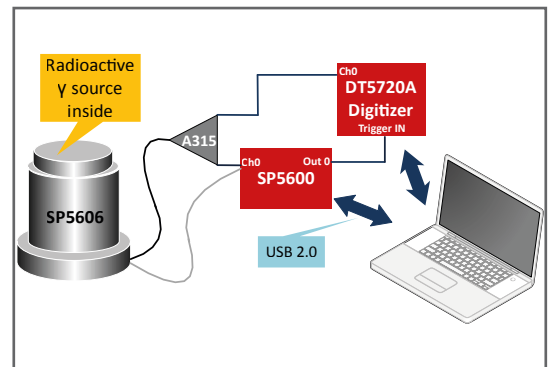

Experimental setup block diagram.

### Results

The student may get acquainted with the presence of radioactivity with a simple preliminary measurement, namely comparing the counting frequency as a function of the discriminator threshold with/without the source. Presuming the source, essentially in contact to the crystal, to be point like with respect to the crystal surface, and assuming its activity is known, the student may estimate for every threshold value the detection efficiency and the signal over noise ratio, building up an efficiency-purity plot. Exemplary results obtained with a  $^{137}\text{Cs}$  source are shown. Moving away the source from the crystal, the law governing the variation of the flux can also be investigated.

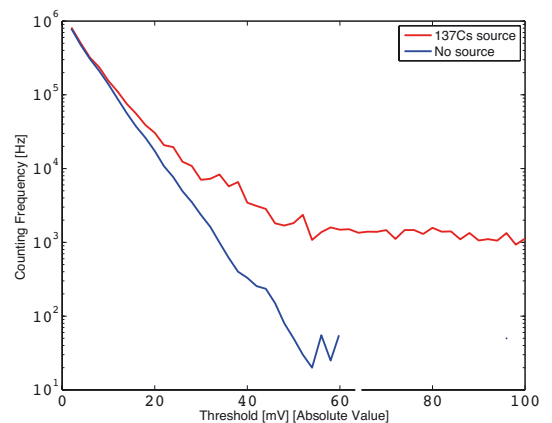

Sensor output frequency as a function of the threshold in mV, with and without  $^{137}\text{Cs}$  source.

This experiment is also possible with the following kits

**B** SP5600AN  
Educational kits  
★ PREMIUM



see  
p. 179

**C** SP5640  
GammaEDU



see  
p. 183

**D** SP5630EN  
Environmental kit



see  
p. 181

**E** SP5630ENP  
Environmental kit Plus



see  
p. 181

**F** SP5600EMU  
Emulation kit



see  
p. 180