

C.1.10 Environmental and Cosmic Radiation

SG6218D



Dedicated kit	
Description	pp.
SP5620CH Cosmic Hunter	182



Requirements

No other tools are needed.

Difficulty	Execution Time	Data Analysis	Radioactive Sources
		YES	NO

Equipment

SP5620CH - Cosmic Hunter

Model	SP5621	SP5622 (x2)
Description	Coincidence Module 	Detection System 

p. 182

p. 187

Purpose of the experiment

To estimate the contribution of environmental radiation during the detection of the cosmic radiation.

See the
Application




Fundamentals

Natural radiation is composed of two components: Environmental (primarily derived from soil, but also water, air and food) and Cosmic radiation. Gamma rays coming from soil will have mean energy lower than 2MeV and will descend from the decay chains of three natural radioactive elements: Potassium [^{40}K], Thorium [^{232}Th] and Uranium [^{238}U]. However, there is a not null probability that a soil-derived gamma interacts with the scintillating tile and deposits 1MeV of energy. This potential could cause confusion or inaccuracies in cosmic rays counting measurement. However, the high threshold of the electronics and the low probability of this phenomenon, when compared to the cosmic ray high detection probability, should typically avoid this inconvenient situation.

Carrying out the experiment

Connect the cable connectors of the two SP5622 to the tile inputs located on the rear panel of the SP5621 module. Power on the SP5621 module and start the acquisition via the front panel START button. When a charged particle crosses the black tile its energy is converted into scintillation light. The photons which are produced are detected by the photosensor and converted into an electrical signal. The number of counts for each scintillator may be viewed via the SP5621 display. Select double scintillators coincidence mode via the related button on the front panel, and then select measurement integration time.


Before beginning acquisition the geometry of the system must be determined. Be sure to keep this geometry for the duration of the experiment. To get acquainted with environmental radiation detected by the Cosmic Hunter place the system over bare soil and record data. Once complete, place the detectors a few meters above bare soil and place an iron or lead shield between the soil and the tiles. Acquire the data again with identical tile geometry and compare the results.

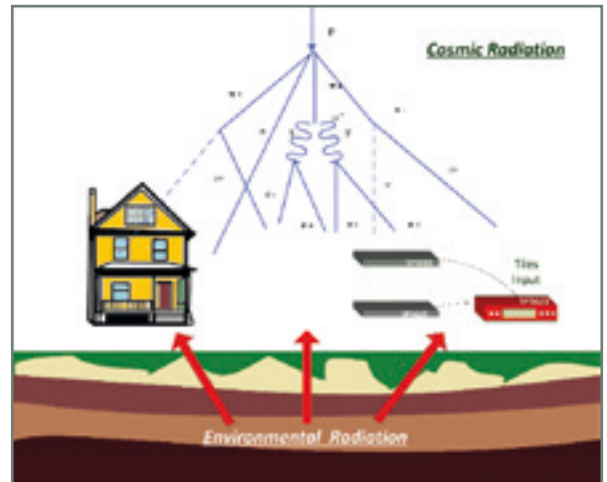

Experimental setup block diagram.

Results

Students may become acquainted with the presence of natural radioactivity by identifying environmental and cosmic contributions via simple comparison of the counting measurements at different heights.

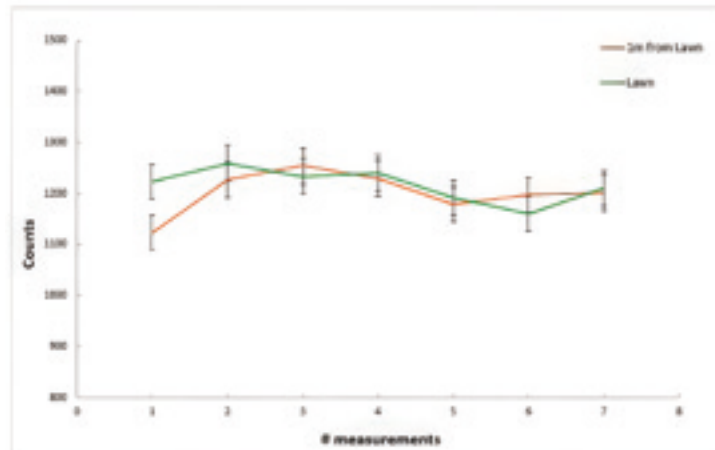

Comparison between the count rates of the measurements acquired on soil and again 1 meter above soil. The comparison plot demonstrates that the parameter settings (Bias Voltage and Threshold) are such that cosmic ray detection is not affected by environmental radiation