



CAEN Educational Handbook

Nuclear and Particle Physics Experiments



<https://edu.caen.it>

Difficulty



Execution Time



SP5630EN - ENVIRONMENTAL KIT



Purpose of the experiment

Get familiar with radon passive measurements by taking care of the proper sample preparation and exposure, as well as of acquiring the spectrum to calculate the Radon concentration.

Fundamentals

Radon is a naturally occurring radioactive gas produced by the breakdown of uranium in soil, rock, and water. Radon can be dangerous since it accumulates inside houses or buildings. Air pressure inside your home is usually lower than the pressure in the soil around the building foundation. Because of this difference in pressure, the

building acts like a vacuum, drawing radon in through foundation cracks and other openings. Additionally, building materials (such as granite and certain concrete products) can give off radon.

It is usually recommended to make screening measurements to have a quick estimate of the highest concentration and take action in case the measurement exceeds 4 pCi/L (or 100 Bq/m³), which is the limit for a non-dangerous exposure to this element.


Requirements

No other tool is needed.

Carrying out the experiment

Detailed instructions for this activity are reported in GD8359 - SP5630EN QuickStartGuide.

Check the expiration date of the charcoal sample; if expired, or if the sample has already been used, bake it for some hours. Acquire the background spectrum with the i-Spector. Before starting the acquisition, make sure that the temperature is stable from the web interface. Once done, expose the sample for about 5-6 days, following the requirements of the closed house, and far away from wall, windows, etc., according to the Radon measurement procedure. Once ready, seal again the sample, wait at least three hours, and acquire the energy spectrum. Take few hours of acquisition with the exposed sample in the same condition of background acquisition.

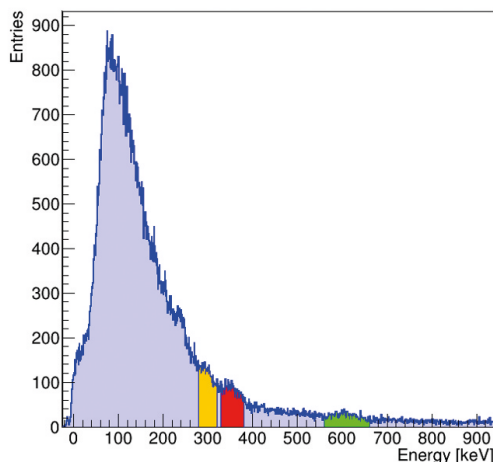

Radon spectrum: counts of the three ROI are summed and combined with the calibration curves. The result is 124 ± 85 Bq/m³. Longer acquisition time will improve the measurement resolution.



Experimental setup.

Results

After the energy calibration and background subtraction, it is possible to select the region of interest (ROI) of the ²¹⁴Pb and ²¹⁴Bi decay products at 295 keV, 352 keV, and 609 keV. The counts can be used together with the calibration curves to calculate the Radon concentration. On request, Excel® spreadsheets examples can be provided to customers.